



Instructions for use

KIMTECH* Wipers for the WETTASK* System

STEP 1

Insert a KIMTECH* Wiper roll into the bucket or canister. Fill out the label with all required information.

STEP 2

Mix your preferred solution for cleaning and/or disinfecting and slowly pour the chemical over the top of the roll.

STEP 3

Thread the centre wiper through the lid and snap shut.



Note:

- Understand any hazards and wear correct Personal Protection Equipment
- Only refill the container with the same wiper code and chemical
- Recommended liquid amount to be added: Bucket 1.9 L/Canister 800ml